

Isohelix
DNA Sampling and Purification


GeneFiX

Saliva DNA/RNA Collection

STORE AT ROOM TEMPERATURE

www.isohelex.com

15°C / 30°C

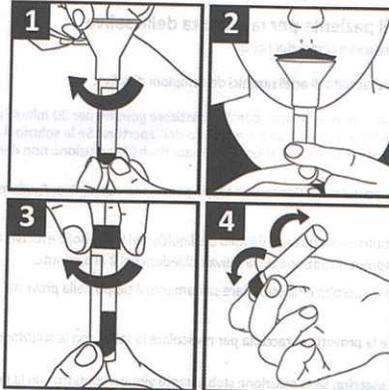

REF GFX-02 / GFXH1-02 / GFXH2-02

Cell Projects Ltd
2 Roebuck Business Park,
Maidstone, Kent, UK, ME17 1AB

RESEARCH USE ONLY

Patient Instructions for Saliva Collection

ENGLISH

Read all instructions prior to collection.

Intended use: For the collection of nucleic acids from saliva samples.

Do not eat, drink, smoke, brush your teeth or chew gum for 30 minutes before use. Check fluid level is at 2ml before opening. If solutions appears cloudy before use, warm tube between hands until solution becomes clear.

1. Remove the cap from the tube. Keep cap for later use. Screw collection funnel on to the tube.
2. Spit into the funnel until the level of the liquid in the tube, not including bubbles, reaches the 4ml line. Delivery of saliva will take between 2 and 5 minutes.
3. Unscrew the collection funnel and replace the tube cap tightly. Discard the funnel.
4. Shake the collection tube several times to mix the saliva with the solution.

Precautions: Do not swallow. If stabilizing solution comes into contact with skins or eyes, wash with plenty of water. Tube cap may be a choking hazard to small children. MSDS available at www.isohelex.com

Contents: Funnel, Tube Pre-Filled with Stabilization Buffer, Cap.